



Delivering Solutions™

◆ Nutritional ◆ Enrichment ◆ Medicated ◆ Special Needs

Product# F7015

Acacia Gum, NF/FCC, 1 kg/Bag

Mixing Directions for Preparing Acacia Gum for Use with Primate Enrichment Devices

1. Add 3 parts powder with 1 part warm water.
2. Mix step 1 by hand or with blender.
3. Acacia gum should be a moist and sticky consistency, but not runny, for easier application.
4. Bio-Serv's flavored PRANG™ can be substituted for the water to provide added variety.